


# MEA-STIM-ADPT-UP-BC

## Pin layout

Adapter for MEAs with 16 additional stimulating electrodes (MEA-200/30-Stim)  
and for MEA1060-UP-BC amplifiers

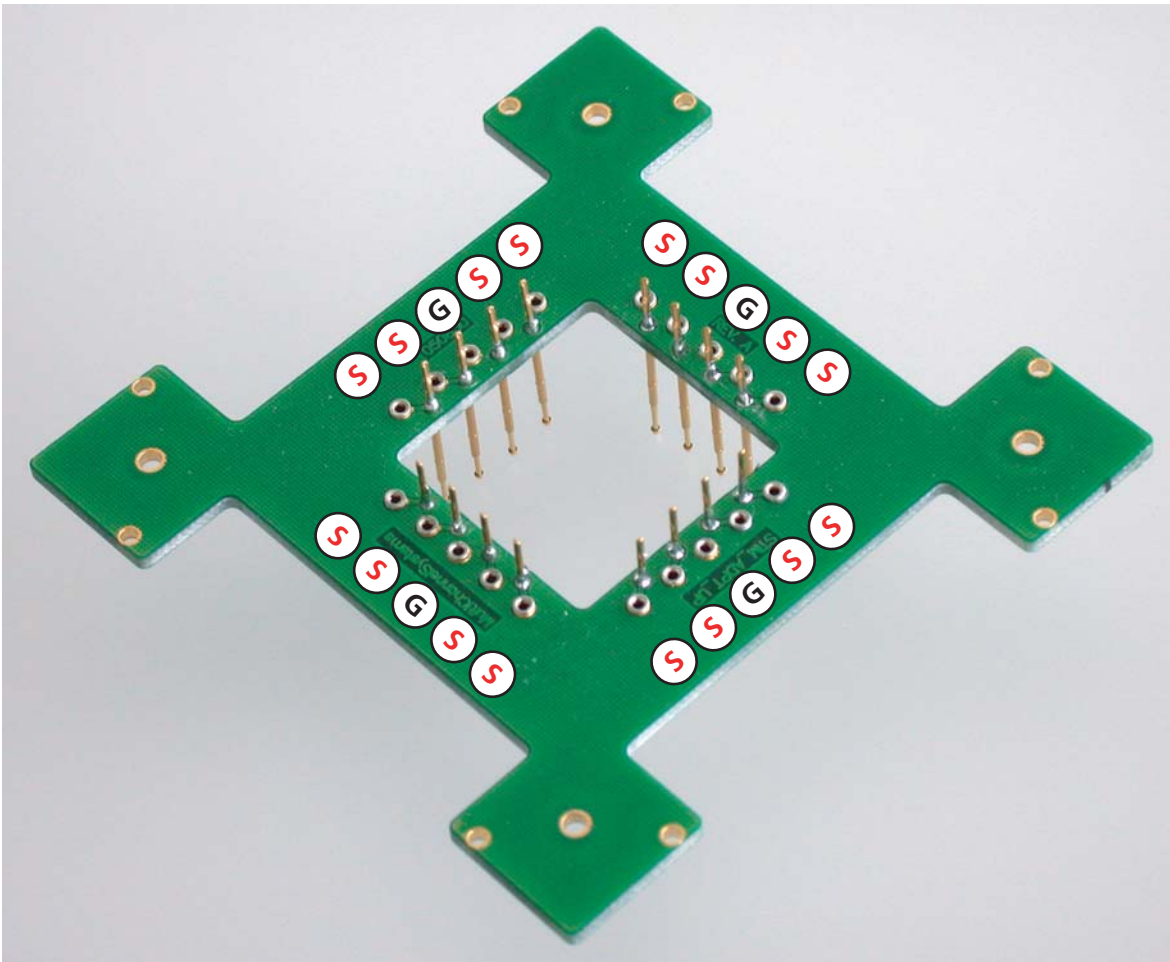

ⓐ = Ground    Ⓢ = Stimulating channel

The adapter is installed on the MEA-1060-UP-BC amplifier as follows. Firstly, the four screws on the amplifier's lid are removed. Then, the adapter is placed onto the four mounting holes, and then tightly secured with the screws. The adapter is symmetrical, that is, the orientation does not matter.