
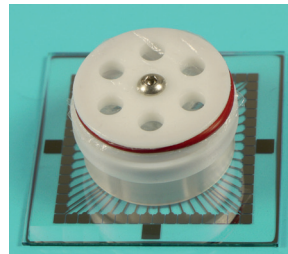


## 6well-CC

### Culture Chamber for 6wellMEAs with triangle or round chamber rings

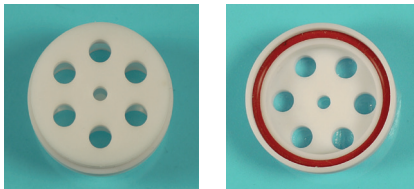

The 6well-CC culture chamber from teflon for mounting on MEA rings with triangle or round chambers is fixed with a socket head screw.

### Application

The culture chamber 6well-CC for 60-6wellMEAs or 256-6wellMEAs is suitable for the 6 well round chamber ring as well as for 6 well triangle chamber ring. To use a 60-6wellMEA or 256-6wellMEA in combination with the 6well-CC as a culture chamber, please insert an O-ring inside the 6well-CC for sealing. Cover the culture chamber with foil and fix the foil with a second O-ring. Then mount the 6well-CC on the triangle or round chamber ring on the MEA. Make sure to place the holes exactly above the chambers of the MEA ring. Fix the culture chamber 6well-CC carefully with a socket head screw on the 6 well ring glued on the MEA.

**Important** The foil to cover the 6well-CC has to be ordered separately!

For example: ALA MEA-MEM-SHEET (Set of 10 3" x 3" sheets of fluorinated ethylene-propylene membranes for sealed MEA culture dishes). Can be ordered at Multi Channel Systems MCS GmbH.

Scope of delivery:	1	6well-CC
	2	O-rings (26 x 2 mm)
	1	Socket head screw (M3 x 12)

### Additionally available from Multi Channel Systems MCS GmbH

6well-CC-Set  
Set of five 6wellCC culture chambers for 6wellMEAs

Scope of delivery:	5	6well-CC
	10	O-Rings
	5	Socket head screws (M3 x 12)
	1	Allen key
	1	Box of membrane sheets (ALA MEA-MEM-SHEET, 10 sheets of 3 x 3")

