


ADPT-FM-32

FlexMEA36 Adapter for ME2100-HS32 and 32-Channel Miniature Preamplifiers MPA32I

Example:
Connecting the FlexMEA36 via
Adapter ADPT-FM-32 to the MPA32I.

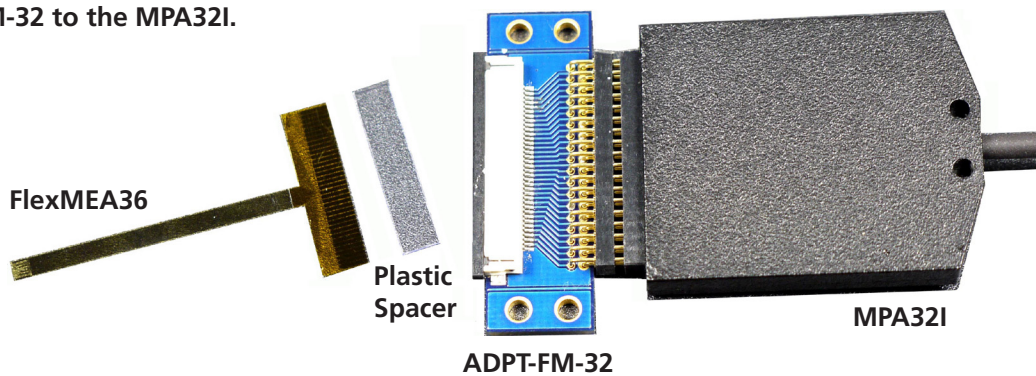

Application

MPA32I: The side with no screws is considered the top side of the MPA32I, please see the foto.

ADPT-FM-32: If the adapter ADPT-FM-32 is oriented as shown on the foto above the FlexMEA36 must be inserted with the contact pads facing downward. Open the white drawer of the adapter.

FlexMEA36: Insert the FlexMEA36 carefully into the drawer of the adapter. Please use the plastic spacer to fix the FlexMEA36 in the drawer of the adapter ADPT-FM-32 to ensure a good contact.

Plastic Spacer: Insert the plastic spacer carefully into the drawer of the adapter. The plastic spacer should be placed on the opposite side of the contact pads of the FlexMEA36, otherwise the connection will be interrupted.

Close the white drawer of the adapter ADPT-FM-32.

If you like to use the FlexMEA36 together with the ME2100-HS32 headstage or with the standard 32-channel miniature preamplifier MPA32I, you need the ADPT-FM-32 adapter to connect the FlexMEA36 to the standard MPA32I. There is no need for an adapter if you use the 32-channel miniature preamplifier MPA32I-Flex, which is specified for FlexMEAs. Please read the MPA32I-Flex manual for more information.