


● **Waste Bottle**  
for Constant Vacuum Pump CVP  
with Pressure Control

● **Instructions for the Replacement**  
of the Waste Bottle of the CVP



**Waste Bottle**

**Top Cover with Tube Connector**

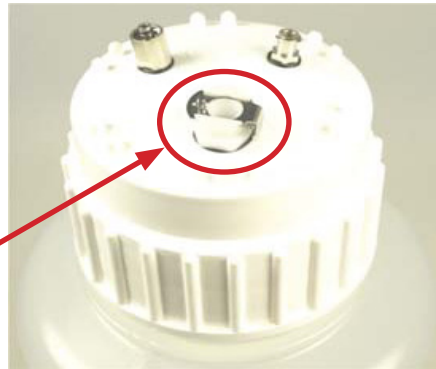

● **Application**

● Please connect the connector of the provided tube to the designated connector on top of the cover of the waste bottle.

Insert the waste bottle into the fitting of the constant vacuum pump. Turn the bottle until the open connectors of the cover are in the back. Connect the tubes which are fixed inside the CVP to the connectors on top of the waste bottle. Connect the right tube to the right connector and the left tube to the left connector.

The capacity of the waste bottle is 5 liter.



**Tube with Connector**