


## CMOS-TH

## Tissue Holder for CMOS-MEA5000-System

### Cleaning

Rinse the tissue holder with distilled water.


Rinse shortly with alcohol for sterilization.

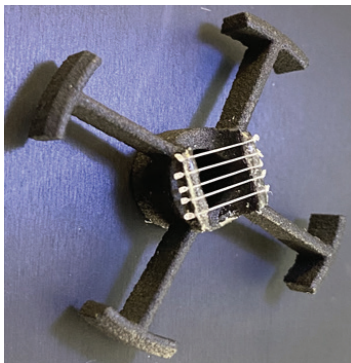

Tissue holder, bottom side with grid

### Scope of delivery

1 CMOS-TH

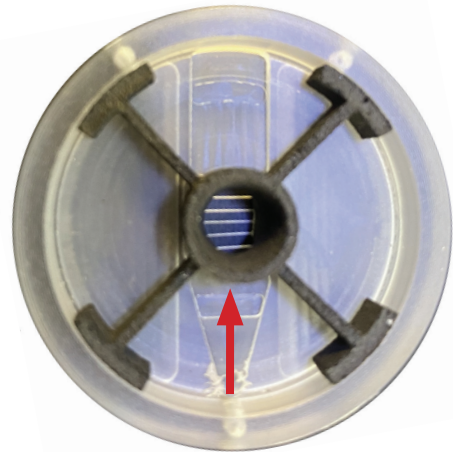

Tissue holder inserted in a CMOS chamber

### Tissue Holder for CMOS-MEA5000-System

Tissue holder for fixing the slice in the CMOS MEA chamber on the electrode field for better recording results.

The CMOS-TH is adapted for the slice chambers on the CMOS chip [Slice Chamber Basic SCB](#) and [Slice Chamber Advanced SCA](#). Both chambers have a flow direction for the perfusion fluid and the tissue holder should be orientated in the same direction.

Please see the arrow in the picture above.

**Important:**  
**Do not autoclave the CMOS-TH tissue holder!**