


Replacement Instructions for the Foil  
of the Modular Setup of the PAL24/96

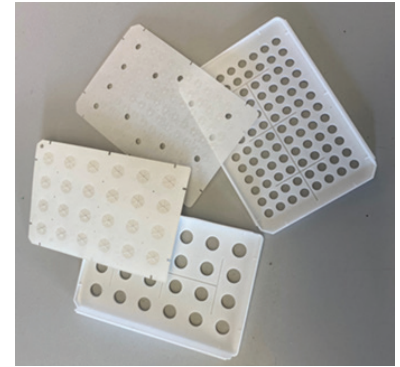


## Replacement Foil for the PAL24/96

Technical Specifications	
Operating Temperature	10 °C to 40 °C (40 °F to 104 °F)
Relative Humidity	20 - 80 %
Foil	
Number of covered wells	24 or 96 wells
Inlet connection to CO2 tube	0.9 mm luer connector (braun sterican cannula)
Dimension (W x D x H)	85.48 mm x 127.76 mm x 0.3 mm
Foil element	Polyurethan
Number of cross cuts for wells	24 o 96

### Mounting Instructions for the Foil of the PAL24/96

1. Remove the old Polyurethan foil from the PAL24/96 lid.
2. Clean the PAL24/96 lid with warm water and a mild detergent. Dry the lid completely.
3. Provide the new foil for the respective lid, PAL24 or PAL96.

The foil has to be fixed at the bottom of the lid.

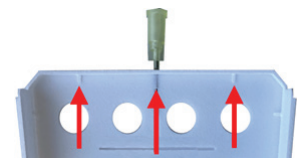

4. Peel off the protective sheet for 2 cm from the foil. Crimp the protective sheet as shown on the picture on the left.
5. Locate the foil precisely that the slots of the foil meet the noses of the lid. Please see the picture above.
6. Subtract the complete protective sheet. All slots of the foil have to meet the noses of the lid. Fix the foil carefully with a weak shim.

