


MEA1060-Up-BC-PGP

Perfusion Ground Plate
for MEA1060-Up-BC Amplifier

Adapter to use perforated MEAs (pMEA)
together with MEA1060-Up-BC Amplifiers

The external diameter of the cannula Perfusion In and Perfusion Out is 0.9 mm. It is recommended to connect flexible hoses with an internal diameter of 0.8 mm that enclose the cannula leakproofed.

Note: Please do not clean the plexiglass of the perfusion ground plate with alcohol! The plexiglass will get refractory. Please use distilled water only.

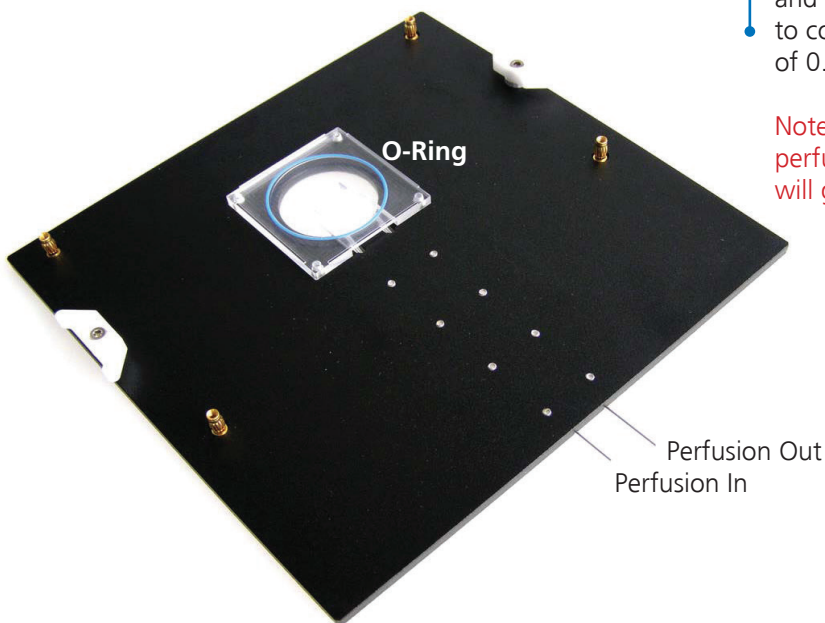

Installation

Replace the standard ground plate of the MEA1060-Up-BC amplifier by the MEA1060-Up-BC perfusion ground plate. Please note that the MEA1060-Up-BC-PGP does **not** have a heating element.

Scope of delivery

- 1 x MEA1060-Up-BC-PGP
- 3 x Silicon sealing rings (PGP-O-Ring 40 x 1, available in units of four rings)
- 1 x Set of screws